


INSTRUCTION MANUAL FOR MAGNISCOPE VIEWER MG-600

Thank you for purchasing the ELMO Magniscope Viewer MG-600.

The MG-600 is designed to magnify glass slides or micro slides in conjunction with either the model EV-4400AF, HV-5000XG, EV-2000AF or HV-3000XG Presenter. (Optional light viewer is needed for the use with EV-2000AF & HV-3000XG)

Direction for use

1. Place the MG-600 on the stage of the Presenter.
2. Turn the camera head of the Presenter to the bottom, fit the camera head to the lens part of MG-600. Attach the close-up lens on the Presenter at this time.
3. Turn on the base light of the Presenter or turn on the light viewer (optional) for EV-2000AF & HV-3000XG.
4. Set the lens of the Presenter to maximum telephoto position.
5. Turn the focusing knob of the MG-600 to maximum far position.
6. Place the prepared slide on the stage of MG-600.
7. Turn the focusing knob of the MG-600 to adjust to the sharpest image.

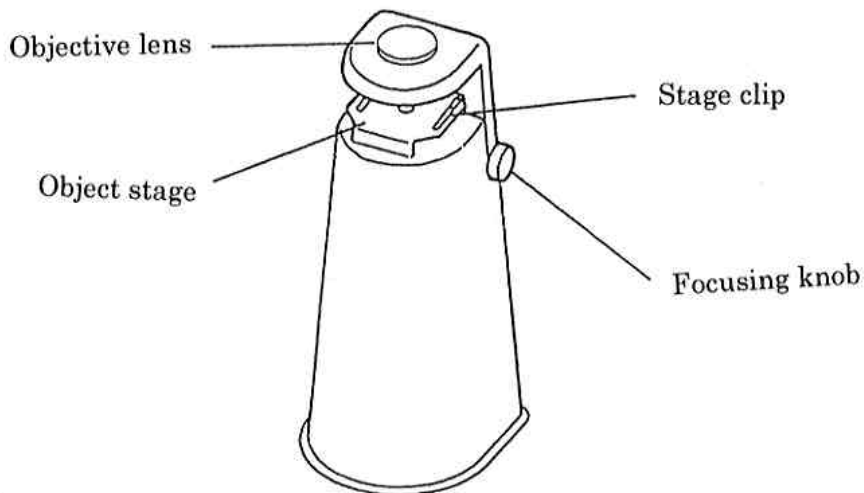

Specifications

Dimensions : H 355mm (14.0") x W 152mm (6.0") x D 157mm (6.2")

Weight : Approx. 900grm (2.0lbs)

Illumination : Utilizes the built-in baselight of the Presenter or the optional light viewer.

Magnification : Approx. 200 times with EV-4400AF

Approx. 130 times with HV-5000XG

Approx. 170 times with EV-2000AF

Approx. 150 times with HV-3000XG (when viewed on 20" monitor)

***ELMO* ELMO CO., LTD.**

6-14, Meizen-cho, Mizuho-ku, Nagoya, 467-8567 Japan E-mail: foreign-div@elmo.co.jp

ELMO Mfg. Corp.

1478 Old Country Road,
Plainview, NY 11803-5034, U.S.A.
Tel. 516-501-1400 Fax. 516-501-0429
E-mail: elmo@elmousa.com
Web: <http://www.elmousa.com>

ELMO Canada Mfg. Corp.

44 West Drive, Brampton,
Ontario, L6T 3T6, Canada
Tel. 905-453-7880 Fax. 905-453-2391
E-mail: info@elmocanada.com
Web: <http://www.elmocanada.com>

ELMO (Europe) G.m.b.H.

Neanderstr 18,
40233 Düsseldorf, Germany
Tel. 0211-376051-53 Fax. 0211-376630
E-mail: elmo@europe@AOL.com
Web: <http://www.elmo.de>

Printed in Japan