




ELMO TRAINING SESSION

USING A MICROSCOPE WITH THE ELMO TT-02U

(Note that this covers all ELMO TT products that do not have a LED Light)
(TT-02 and TT-02U only)


Teachers who use a Microscope will find that the ELMO TT-02 or TT-02U document camera when attached LCD projector is an amazing tool. Imagine that finally you can show the entire class an image from the Microscope at the same time. No other accessories are needed!

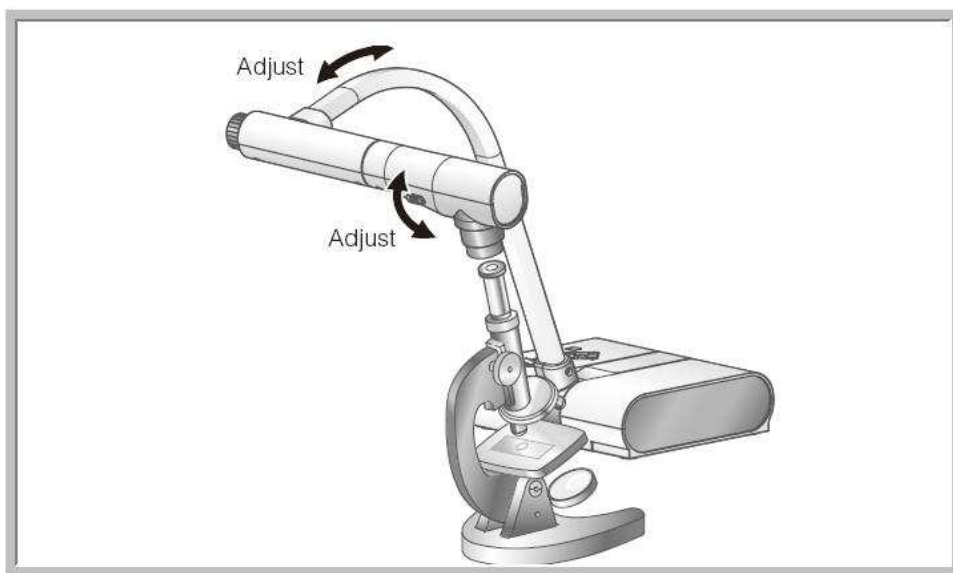


Follow these easy steps

- Move the neck and camera of the TT-02 or TT-02U so that the camera is looking down the eye piece of the microscope. Make sure that the TT-02 or TT-02U is Not Touching the eyepiece of the Microscope at this time due to the fact that the TT-02 and TT-02U lens moves out when setting the Microscope Macro. You do not want to put any pressure on the document camera lens.


- Click on the Menu button

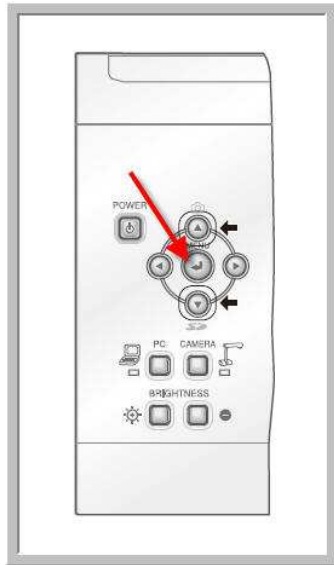

Then click on the top or bottom arrow until you see the word "Microscope"

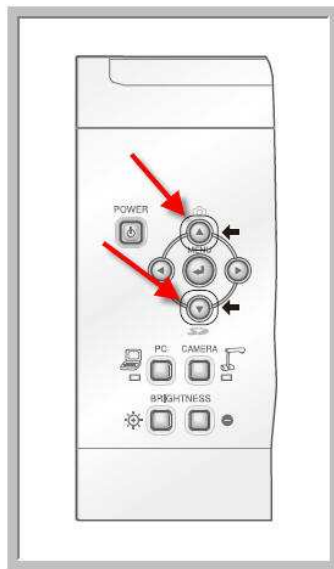

Click on the right button (see below) and highlight “SET” and hit the Menu button

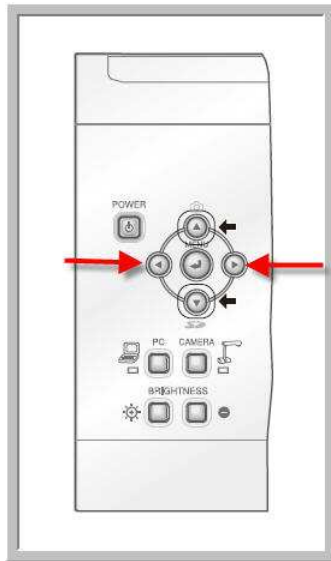

- You have just set the “Macro” mode to the ELMO TT-02 or TT-02U. Now place the Elmo document camera lens on top of the microscope eyepiece. Now you have to fine tune the image using the Microscope to Zoom and position the image. Note that you might have to recenter the camera lens of the ELMO document camera.
- Finally, if you have an ELMO Document camera with a USB (TT-02U) you can store the image shown on the microscope if your are attached to a Windows computer via USB. ELMO Image mate software has to be previously loaded on the computer.
- Remember to reset the Elmo TT-02 or ELMO TT-02U to Microscope “Off” mode by using the menu button and the microscope setting when you want to view documents or anything other than from the microscope.

If you have any questions please email me at agraziano@elmousa.com.